DECLARACIÓN DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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*This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.*

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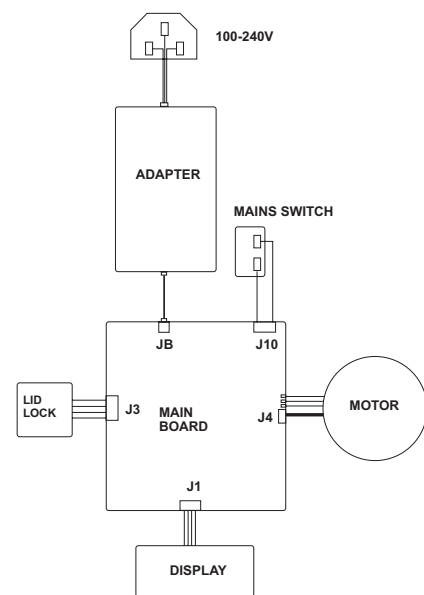
**ANEXO I: WIRING DIAGRAM FOR CENTRIFUGE**

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

**INSTRUCTIONS ON ENVIRONMENT PROTECTION**

Do not dispose of this product in the usual household garbage at the end of it's life cycle; hand it over at a collection point for the recycling of electrical and electronic appliances. The symbol on the product, the instructions for use or the packing will inform about the methods for disposal.

The materials are recyclable as mentioned in it's marking. By recycling, material recycling or other forms or re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

**7. TROUBLESHOOTING**

ERROR	CAUSE	SOLUTION
No display	No main power connection Power failure	Plug in mains cable on both sides Check the mains fuse
Lid can't be opened	Power failure. Rotor is still spinning. Lid locks failure	Emergency lid release (see sec.4.7) Stop centrifuge run Contact service
Centrifuge shakes during acceleration	Rotor not loaded symmetrically	Stop centrifuge run and load centrifuge symmetrically.
Er=01	Open lock failure	Contact service

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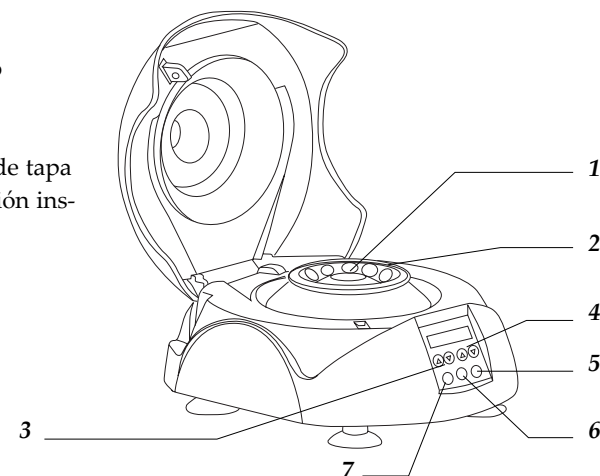
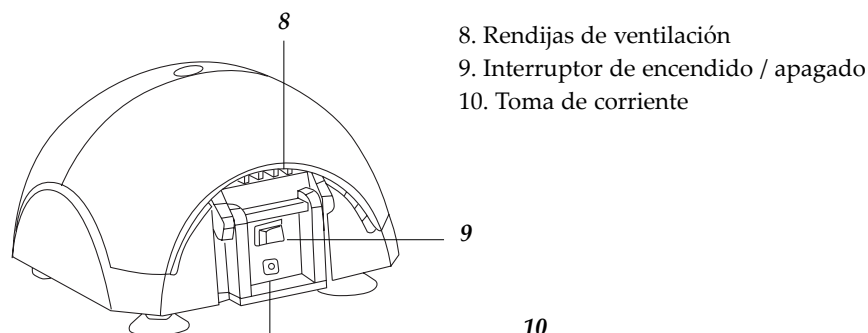
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

### 6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ Even slight scratches and tears of the rotor can lead to serious internal material damage.
- ◆ Avoid damage caused by aggressive chemicals, including among others: strong and weak alkali, strong acid. In the event of contamination caused by aggressive agents, the rotor must be cleaned immediately using a neutral cleaning liquid.
- ◆ The equipment has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

#### Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the equipment and produce an early ageing of the equipment.
- ◆ The rotor and the outside of centrifuge should be cleaned regularly with a fluff-free cloth dampened with soaped water that does not contain abrasives. Disconnect the centrifuge from the mains supply, remove the rotor and clean it separately. Only neutral agents may be used for cleaning.
- ◆ Please check the rotor and especially the rotor bores regularly for deposits or damage, then reinsert the rotor and tighten the rotor nut.

to end the run before the time finish. The remaining run time appears in the display.

After the run, the lid lock of centrifuge opens automatically.

The time setting and the rotational speed may be changed during the run. The remaining run time appears in the display.

#### 5. Short-spin centrifugation

- ◆ Turn on the mains switch, pressing "OPEN" to open the lid, load the rotor symmetrically. Fasten the rotor lid and close the centrifuge lid.
- ◆ Pressing "SHORT-SPIN". Short-spin centrifugation is possible for as long as this key is held down. The maximum rpm is 13400.

#### 6. Switching to the rcf display

- ◆ Pressing speed's ▲▼ simultaneously, the display switches from rpm to rcf. Also pressing speed's ▲▼ simultaneously again, the display switches from rcf to rpm.

You can calculate with the following formula :

$$RCF = 1118 \times 10^{-5} \times n^2 \times r_{max}$$

n: rotational speed in 1/min

rmax= 6 cm, max. centrifuging radius in cm

#### 7. Open the centrifuge in the event of a power failure

- ◆ Disconnect the centrifuge from the mains supply. Wait until the rotor has come to a standstill, then left up the device, remove the lock pole to right, the lid lock will be opened.

**NOTE: If you are not using the equipment for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).**

#### Security

- ◆ The centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ You should put the equipment in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ◆ Do not place the centrifuge near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ◆ When you are not using the equipment for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).

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#### *Setting up*

1. Control elements
2. Rotor insertion and removal
3. Loading the rotor
4. Centrifugation with a preset time
5. Short-spin centrifugation
6. Switching to the rcf display
7. Open the centrifuge in the event of a power failure

##### *1. Control elements*

START/STOP: for start or stop centrifugation

SHORT-SPIN: for reduced-time centrifugation

OPEN: to open the lid

▲▼: for setting the time and speed

##### *2. Rotor insertion and removal*

◆ Fit the rotor onto the motor shaft, followed by the rotor nut. Tighten the rotor nut by turning clockwise. To release the rotor, turn the rotor nut counter clockwise.

**IMPORTANT** Before each start, check that the rotor is firmly tightened!

◆ The rotor and the rotor lid must always be securely fastened. If the centrifuge makes unusual noises when started, the rotor or rotor lid is not fastened correctly. Switch the device off immediately by pressing "START/STOP" key.

**IMPORTANT** Do not use damaged rotors!

**IMPORTANT** Do not move the centrifuge during the run!

##### *3. Loading the rotor*

◆ Rotor must always be loaded symmetrically. Minimize differences in weight between the filled sample tubes-taring with a scale is recommended. This will reduce wear on the drive and cut running noise.

◆ A liquid density of 1.2g/ml must not be exceeded at the max. speed.

◆ Damaged tubes can not be centrifuged. This is because broken tubes can, in addition to sample loss, result in further damage to the centrifuge.

◆ Close the test tube lids before centrifuging. Open lids can be ripped off during centrifuging and damage the centrifuge.

##### *4. Centrifugation with a preset time*

◆ Turn on the mains switch, pressing "OPEN" to open the lid, load the rotor symmetrically. Fasten the rotor lid and close the centrifuge lid.

◆ Pressing Time's ▲▼ change the run time, the time can be pre-selected between 20 seconds and 99 minutes, and pressing speed's qp to change the run speed, the max. speed is 13,400rpm.

◆ Pressing the first "START/STOP" to start the run. Pressing the second "START / STOP"

## 5. INSTALLATION / SETTING UP

### Preliminary inspection

Unwrap the centrifuge, take off the involving plastic and take off the polispán protection in which it comes fitted. Take off all the protective items and, without connecting the equipment to the net, make sure that it does not present any damage because of the shipment. In case the equipment presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Rotor
- Rotor nut
- Rotor lid
- Adapter
- Operations manual
- Warranty certificate

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

### Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

### PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT

- ◆ Please put the centrifuge on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Make sure that the ventilation slits are not blocked.
- ◆ Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ Avoid inflammable or toxic substances in the working area.
- ◆ Insert the column connector of the adapter to power connector of the device, and insert another connector of the adapter to mains power supply.
- ◆ Power on the main switch. The centrifuge is ready to operate when the display becomes visible. Place the rotor onto the rotor axle and tighten using the rotor nut.

**IMPORTANT: Before starting up the centrifuge for the first time, make sure the rotor nut is securely fastened.**

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz  $\pm$  10%.

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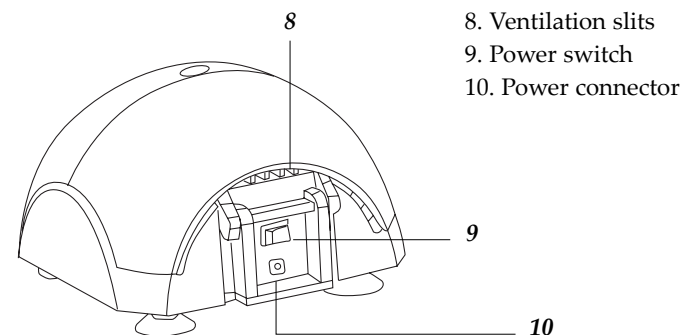
### 4. THEORETICAL FOUNDATIONS

Centrifugation is a separation technique based on the movement of the particles that are displaced to the distal end of the rotation axis depending on their shape and mass.

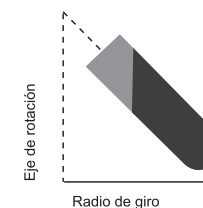
The Relative Centrifugal Force (R.C.F.) is the required force to separate particles. It is given as some value times that of gravity (g).

$$R.C.F. = 1.118 \cdot 10^{-5} \cdot r \cdot n^2$$

$1,118 \cdot 10^{-5}$  is a constant.

$r$  = radius, the horizontal distance in centimetres from the rotation axis to the bottom of the sample tube.

$n$  = rotation speed expressed in revolutions per minute (r.p.m.).

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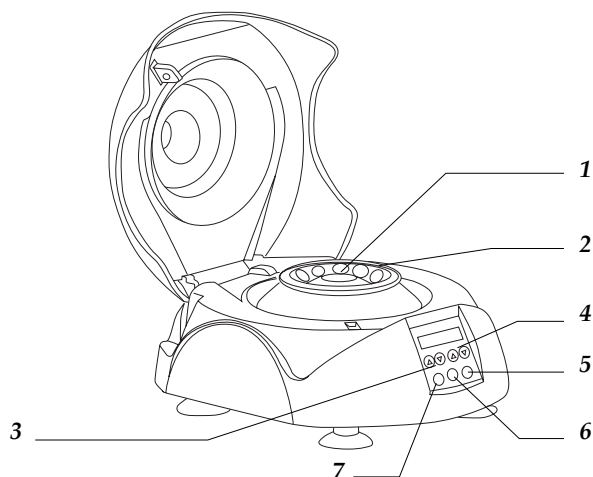
**1. USES OF THE INSTRUMENT**

This centrifuge is designed for workstations in training and research laboratories in the fields of bioscience, medicine and chemistry. Twelve 1.5ml or 2.0ml micro test tubes can be centrifuged simultaneously in a 45° fixed-angle rotor.

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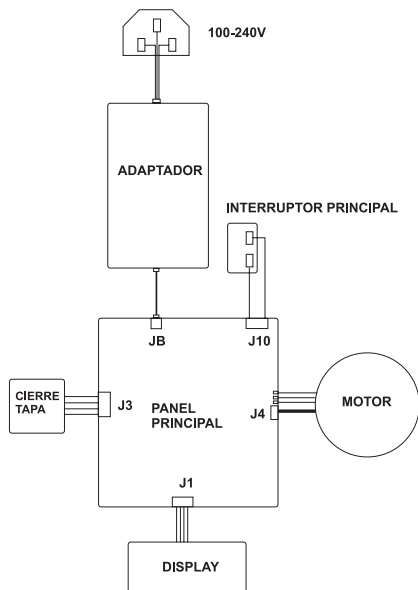
Thank you for choosing this equipment. We sincerely wish that you enjoy your centrifuge Nahita 2507/100. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the centrifuge Nahita 2507/100, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the equipment.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es) or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the equipment warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment